



## **Judd Rice Laboratory**

*at the USC/Norris Comprehensive Cancer Center*

[www.histonecode.com](http://www.histonecode.com)

### **Annealing Two Single-Stranded Oligos**

Revised by Siddhartha / Trun Lab

August 2003

1. Heat equal amounts of both oligos to 94°C for 5 min.
2. Cool the oligos 1° per 100 seconds (or as slow as the PCR machine will allow).
3. It should take 3.5 hrs to cool.
4. Check the annealing reaction by capillary electrophoresis.